

# HOT FOOT® DOUBLE-LOCK LIGHTER INSTRUCTIONS

## SPECIFICATIONS

Piezoelectric ignition

Double-lock design fulfills U.S. Government Safety Regulations

Butane refill

## OPERATION

1. Slide the ON/OFF switch on the side of the unit to ON.
2. Slide the small black slide switch on top of the unit to the left to reduce the size of the flame; slide it to the right to increase the size of the flame.
3. Press and continue to HOLD DOWN the round black button on top of the unit and pull the trigger.
4. Release the trigger to stop lighting.
5. To refill the unit always hold the butane fuel can upside down. Press the butane can tip vertically downward into the center of the refill valve (found on the end of the lighter) and apply firm pressure.

**CAUTION:** DO NOT refill near an open flame.

**Butane Refill Can** - order butane can (provides for thousands of lights), 2.5 ounces

**Control Company Cat. No. 3402**

For more information, please contact us:

[ExpotechUSA](#)  
[10700 Rockley Road](#)  
[Houston, Texas 77099](#)  
[USA](#)

[281-496-0900 \[voice\]](#)

[281-496-0400 \[fax\]](#)

E-mail: [sales@expotechusa.com](mailto:sales@expotechusa.com)

Website: [www.ExpotechUSA.com](http://www.ExpotechUSA.com)