

TRACEABLE® CREDIT CARD TIMER INSTRUCTIONS

SPECIFICATIONS

Display:	3/8" high LCD
Range:	99 minutes, 59 seconds
Resolution:	1 second
Accuracy:	0.01%
Alarm:	Sounds for 1 minute when zero is reached
Features:	Memory to recall last set time Time Out Count up stopwatch Magnet/VELCRO™ to "stick" anywhere Two-year replaceable battery
Size:	3 3/8" x 2 1/8" x 3/16"
Weight:	0.7 ounce

CONTROL BUTTONS

1. Start
2. Stop/Reset
3. Display Input (numbers 0-9)

COUNT UP MODE

1. Press the START button to start timer.
2. Press STOP once to stop or pause timer.
3. Press START again to continue.
4. Press STOP button twice to clear the display.

COUNTDOWN MODE

1. Using the Display Input buttons, enter the desired countdown time.
Example: to set the count down time for 10 minutes, 30 seconds press 1,0,3,0.

2. Press START button to start timer.
3. Press STOP once to stop or pause timing.
4. Press START again to continue.
5. When alarm sounds, press any button to turn off alarm and recall memory.
6. Press STOP button twice to clear memory.

NOTE: Alarm will sound for 1 minute then automatically recalls memory.

NOTE: The credit card timer can be placed anywhere using enclosed VELCRO™/Magnet. Simply remove backing and attach to center back of timer.

ALL OPERATIONAL DIFFICULTIES

If timer does not function properly for any reason, replace the battery with a new high-quality battery (see "Battery Replacement" section). Low battery power can occasionally cause any number of "apparent" operational difficulties. Replacing the battery with a new fresh battery will solve most difficulties.

BATTERY REPLACEMENT

A faint display, incorrect display or no display are all indicators that the battery must be replaced. To replace the battery, push left cover lock while pulling cover out with the tip of a pen or a paper clip. Remove the exhausted battery and place with a new battery. Replace the back of the unit.

WARRANTY, SERVICE, OR RECALIBRATION

For more information, please contact us:

[ExpotechUSA](#)
[10700 Rockley Road](#)
[Houston, Texas 77099](#)
[USA](#)

[281-496-0900 \[voice\]](#)

[281-496-0400 \[fax\]](#)

E-mail: sales@expotechusa.com

Website: www.ExpotechUSA.com