

TRACEABLE® ROBO-THERMOMETER™ INSTRUCTIONS

SPECIFICATIONS:

Display: 1/4" LCD

Range: -58.0 to 536.0 °F
-50.0 to 280.0 °C

Resolution: 0.1° above -20.0° and
below 200.0°

1.0° below -20.0° and
above 200.0°

Accuracy: **Cat. No. 4149**
±1°C -20 to 100 °C
Cat. No. 4349
±0.4°C -20.0 to 100.0 °C

Sampling

Rate: 1 second

Features: Protective sleeve with pocket clip,
8" stainless steel probe, 180
degree rotating digital display,
extra-wide temperature range.

CONTROLS/INDICATORS

°F/°C button and ON/OFF button.

OPERATION

1. Press the "ON/OFF" button to turn the thermometer on.
2. Press the "°F/°C" button to toggle between displaying the temperature in Fahrenheit and in Celsius.

3. Remove the thermometer from the protective plastic sleeve and insert the probe into the material to be measured.

4. The head of the thermometer rotates 180 degrees in order to read the display at various angles.

5. Press the ON/OFF button to turn the unit off.

ALL OPERATIONAL DIFFICULTIES

If this unit does not function properly for any reason, replace the battery with a new high-quality battery (see "Battery Replacement" section). Low battery power can occasionally cause any number of "apparent" operational difficulties. Replacing the battery with a new fresh battery will solve most difficulties.

BATTERY REPLACEMENT

Erratic readings, a faint display, or no display are all indicators that the battery must be replaced.

Remove the battery cap by turning 1/4 turn counter clockwise. Replace exhausted battery with a new 1.5 volt battery. Make certain that positive (+) side is face up. Replace battery cap. Do not over-tighten the cap. Replacement battery Cat. No. 1039.

WARRANTY, SERVICE, OR RECALIBRATION

For more information, please contact us:

[ExpotechUSA](#)
[10700 Rockley Road](#)
[Houston, Texas 77099](#)
[USA](#)

[281-496-0900 \[voice\]](#)

[281-496-0400 \[fax\]](#)

E-mail: sales@expotechusa.com

Website: www.ExpotechUSA.com