


INTERIOR RELAY LOCATIONS

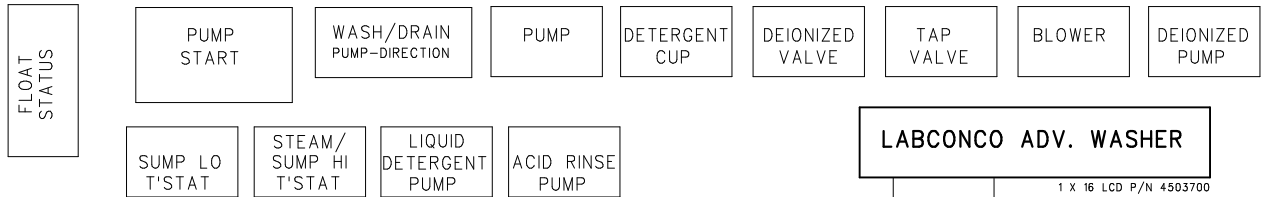


LABCONCO


WIRING DIAGRAM
 ADVANCED WASHER 230V 50/60HZ
 P/N 4534600
 REV. C

WARNING

DISCONNECT FROM POWER SOURCE
 PRIOR TO SERVICING UNIT

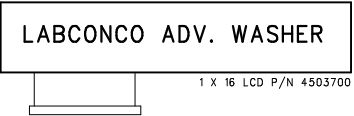



NOTE: PCA FOR REFERENCE ONLY

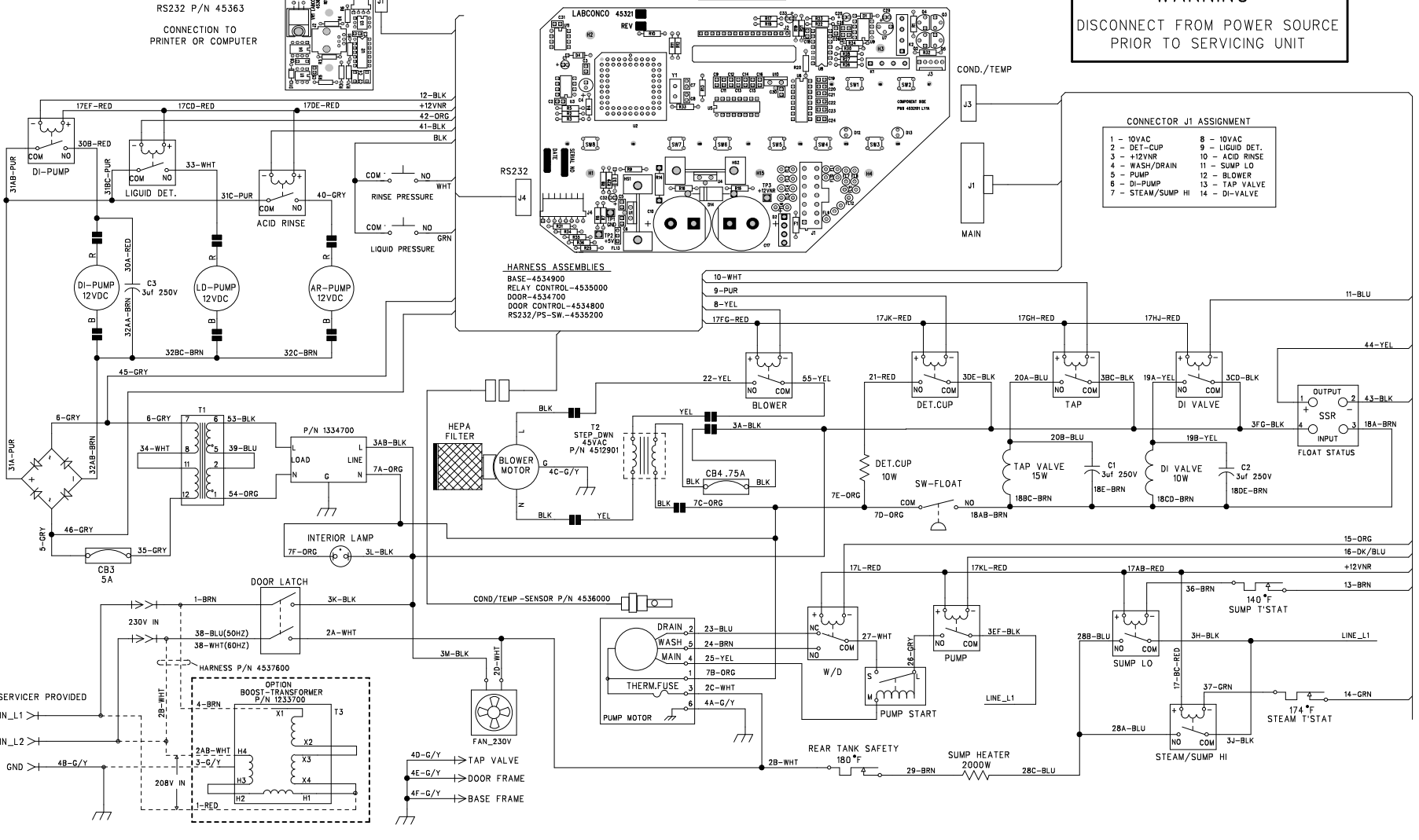

CONNECTOR J1 ASSIGNMENT

1 - 10VAC	8 - 10VAC
2 - DET-CUP	9 - LIQUID DET.
3 - +12VNR	10 - ACID RINSE
4 - WASH/DRAIN	11 - SUMP LO
5 - PUMP	12 - BLOWER
6 - DI-PUMP	13 - TAP VALVE
7 - STEAM/SUMP HI	14 - DI-VALVE

HARNES ASSEMBLIES
 BASE-4534900
 RELAY CONTROL-4535000
 DOOR-4534700
 DOOR CONTROL-4534800
 RS232/PS-SW.-4535200

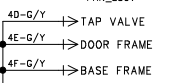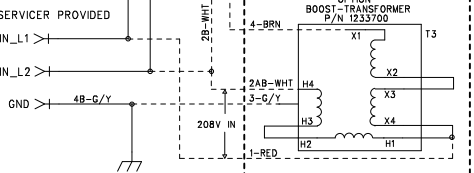

For more information, please contact us:

ExpotechUSA
10700 Rockley Road
Houston, Texas 77099
USA

281-496-0900 [voice]

281-496-0400 [fax]

E-mail: sales@expotechusa.com

Website: www.ExpotechUSA.com