

INSTRUCTIONS:
CHROMALOX 2104-R0000
LOWER/UPPER TEMPERATURE LIMIT SETTING

1. Push “RESET” button for 3+ seconds until “LOCK” appears in upper screen.
2. Then push “UP” or “DOWN” arrow button until “458” appears in the lower screen.
3. Push “RESET” and at the same time, push the “UP” button momentarily until the “INPt” appears in the upper screen and “PAGE” appears in the lower screen.
4. Then push the “RESET” button momentarily “SPLL” – until “SPLL” appears in the upper screen. Then push “UP” or “DOWN” button until 32 appears in the lower screen. This is the lower temperature limit to which the controller can be set for normal operation.
5. Push the “RESET” button again until the “SPUL” appears in upper screen. Then push the “UP” or “DOWN” button until “425” appears in the lower screen. This will be the high temperature limit to which the oven can be set.
6. With these parameters set push the “RESET” button for 3+ seconds and the controller will go back to normal operation with the “SET-POINT” temperature in the lower screen and the measured temperature in the upper screen.

For more information, please contact us:

[ExpotechUSA](#)
[10700 Rockley Road](#)
[Houston, Texas 77099](#)
[USA](#)

[281-496-0900 \[voice\]](#)

[281-496-0400 \[fax\]](#)

E-mail: sales@expotechusa.com

Website: www.ExpotechUSA.com