

NEW

- Compatible with most electrode types
- One-hand operation
- Twistable display

IP 65 CE 3 Years Warranty



pH Meters

VARIO pH

You notice it immediately: apart from its ergonomic form the new VARIO has no keys, but has an innovative touch screen instead. In this way all functions can be called up and set – a simple touch is enough.



VARIO pH

Measuring in no time at all

Just a fingertip touch on the display – and VARIO is ready for use. Immersion in the solution starts the measurement automatically. The stable measurement can be read from the large display together with the temperature and can also be “frozen”. The memory has space for up to 50 measured values for later evaluation.

If the VARIO is not being used for pH measurements then it can be used as a laboratory clock or timer.

Light, handy, robust – it finds a place in every laboratory coat without dripping or leaving nasty stains, as it can be stored without KCl.

The VARIO can operate continuously for 1500 hours on one easily replaceable AA (1.5 V) battery. In addition, the VARIO can operate on line power.

VARIO comes standard with an intelligent glass electrode that is protected by a plastic casing. The conical protective cap does not require KCl which prevents the electrode from dripping and protects it from drying out.

The VARIO can do even more.

The adapter contained in the VARIO Set makes the instrument compatible with commercially available precision electrodes. The VARIO measures as accurately and reliably as a handheld instrument.

You can twist and turn as much as you want, the VARIO is an essential aid whenever speed is required in the laboratory or in production.

Technical Data

	VARIO pH
pH range	-2.00 ... 16.00
pH accuracy	±0.01 pH
Temperature range	-5.0 ... 100.0°C (23 ... 212 °F)
Automatic buffer recognition	TEC/NIST
Calibration points	3 (MultiCal®)

Ordering Information

VARIO	Order No.
VARIO Set	VARIO in the mobile case set, incl. short electrode with built-in temperature sensor and technical buffer 4 and 7 2V00-001V
Other electrodes see brochure “Product Details”	

For more information, please contact us:

[ExpotechUSA](#)
[10700 Rockley Road](#)
[Houston, Texas 77099](#)
[USA](#)

[281-496-0900 \[voice\]](#)

[281-496-0400 \[fax\]](#)

E-mail: sales@expotechusa.com

Website: www.ExpotechUSA.com